



Operating Instructions

Cryofab Pump-out Operator, 1"

To attach, remove plastic cap on pump-out, and push operator over the pump-out body. Push the valve stem down until it touches the plug. Screw clockwise until stem is down, then tighten quick coupling on top. Attach vacuum hose and evacuate the line only. When vacuum pressure in line is down, open the pump-out valve by holding the operator body with one hand while pulling up on the stem with the other.

After evacuation is complete, reseal the pump-out by pushing the stem all the way down. Remove vacuum hose, and then disengage the stem by turning counterclockwise. Loosen quick coupling to make sure stem is free and remove the operator. Replace plastic cap.

Call for information about parts, accessories, and repair services.

1.800.426.2186